

Earthquake Security for your Kuhner shakers



Earthquake Security precautions

Earthquakes can be common occurrences in some countries and Kuhner has taken precautions for the security of your Kuhner incubator shakers. Our machines are designed with built-in stacking pins that shall never detach unexpectedly.


Strong feet

Larger surface area with pre-drilled hole for anchoring to the ground. Reinforced for stability and recommended on all our ISF1-Z triple stacks.



Extra security straps

Strap to walls or floors for added protection

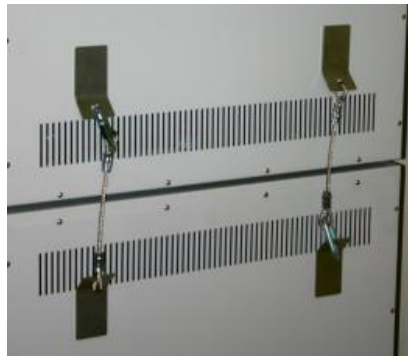

Security rail for water reservoir

Retrofittable aluminum plate with railing screws directly onto the top of our incubator shakers, preventing carboys from being pulled off.



Trust Kuhner for thoughtful engineering, customized projects, easy retrofitting options, and peace of mind for your lab

References: Quakehold Industrial